DECON-AHOL®

WFI Formula/Formule WFI/Fórmula de WFI

For Use in Clean Rooms and Controlled Areas/Pour usage dans les salles blanches et les zones contrôlées/Para uso en salas blancas y áreas controladas
Hard Surface Disinfectant/Désinfectant de surfaces dures/Desinfectante para superficies duras

Active Ingredient/Ingrédients actifs/Ingredientes activos:
Isopropyl Alcohol/Alcool isopropylique/Alcohol isopropílico
CAS#/No CAS/CAS n.°: 67-63-0.....70.0%

Other Ingredient/Autres ingrédients/Otros ingredientes:
*Water/Eau/Agua.....30.0%

Total:.....100.0%

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Veltec Associates, Inc. 15 Lee Blvd. Malvern, PA 19355-1234 USA
1-888-4-STERILE

In Canada, distributed by/Au Canada, distribuído por/

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Canada Clean Room, 200 Terrence Matthews. Kanata, ON. K2M 2C6
Inquiries/Renseignements/ Informes: (888) 595-8070

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USE: Disinfects and sanitizes clean rooms and controlled areas such as those used in healthcare institutions, biopharmaceutical, pharmaceutical, medical device and diagnostic manufacturing facilities. Use on hard non-porous inanimate surfaces in aseptic filling and gowning rooms, general manufacturing areas, or on machinery, tables, counters, laminar flow benches, floors, walls, carts, shelves made of hard non-porous materials, such as glass, plastic, vinyl, chrome and stainless steel. It is compatible with most hard surface materials.

DISINFECTION: Use this product while wearing personal protection. Wear protective eyewear such as goggles or face shield, protective clothing and chemically resistant gloves. Use with adequate ventilation and spray away from eyes and face. **DECON-AHOL** is ready to use. Do not dilute. Pre-clean surface or item of heavy soil or gross filth before application. Hold can upright 6-8 inches from surface. Thoroughly wet surface with **DECON-AHOL** and allow to remain wet for a minimum of 10 minutes. Allow surface to air dry, or after 10 minutes wipe dry with sterilized cloth, wiper, if needed.

DECON-AHOL is bactericidal. The results of AOAC Efficacy Tests show that **DECON-AHOL** is an effective hard surface disinfectant against *Pseudomonas aeruginosa* (ATCC#15442), *Staphylococcus aureus*, (ATCC#6538) and *Salmonella enterica* (ATCC#10708) in the presence of fetal bovine serum organic load in 10 minutes.

SANITIZATION: Use 5 minute contact time on hard, non-porous surfaces (made from: stainless steel, glass, Corian™, laminate flooring, aluminum, pvc, mild steel, polypropylene), in isolators, clean rooms, sterile and controlled areas, gowning rooms and on process lines, gloves, work surfaces, tools and equipment.

DECON-AHOL is an effective sanitizer on inanimate, non-food contact surfaces against *Staphylococcus aureus* (ATCC#6538), and *Klebsiella pneumoniae* (ATCC#4352) in the presence of the fetal bovine serum organic load in 5 minutes.

Storage and Disposal

Do not contaminate water, food, or feed by storage or disposal.

Storage: Store in the original container in a dry cool place below 48.8°C (120°F). Absorb spills with inert dry absorbent or sand. Absorb spills with inert material (e.g., dry sand or earth), then place in a chemical waste container.

Disposal: Product and product wastes are toxic. Follow Federal/Provincial/State regulations and Local/Municipal ordinances when disposing of this product. Improper disposal of excess product, spray mixture or rinseate is a violation of Federal/Provincial/State Laws. If these wastes cannot be disposed of by use according to label instructions, contact your Federal/Provincial/State or Local/Municipal environmental control agency for guidance. Contaminated absorbent in drums must be disposed of via an authorized waste disposal contractor. Never place unused product down any indoor or outdoor drain.

Container Disposal: Non-refillable container. Do not reuse this container to hold materials other than pesticides or diluted pesticides (rinstate). Triple rinse container promptly after emptying. Triple rinse as follows: Fill container 1/4 full with water and recap. Shake 10 seconds. Follow Pesticide instructions for rinstate disposal. Drain for 10 seconds after flow begins to drip. Repeat two more times. Offer for recycling where available or puncture and dispose of in a sanitary landfill, or by other procedures approved by state and local authorities

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